

OptiPrep™ Application Sheets

C47 Isolation of dendritic cells from tissues and cultured cells by sedimentation on to a density barrier

- ◆ **OptiPrep™ is a 60% (w/v) solution of iodixanol in water, density = 1.32 g/ml**
- ◆ For links to other relevant files click on the double blue arrow in the following text

1. Background

Since dendritic cells (DC) were recognized as playing an important role in the induction of cell-mediated responses [1], there has been a rapid growth in research into the function of these cells and methods for their purification. Gradients of either albumin or metrizamide, although providing an effective enrichment of DC, tended to cause some functional alteration of the cells (see ref 2 for details). However, because cells are more tolerant of Nycodenz®, this iodinated density gradient medium rapidly became established as the medium of choice for DC cell purification from peripheral blood and from lymphoid tissues.

More recently methods have been developed for isolation of DC from lymphoid and other tissues by flotation in OptiPrep™ gradients. However in the protocol described in this OptiPrep™ Application Sheet the more traditional method of layering on to a density barrier is described; its merit being that comparison with older Nycodenz® preparations may be easier. This separation protocol has been used for isolation of DCs from lymph [3], bone marrow [4], bone marrow cell cultures [5], spleen [6], thymus [6] and lung [7].

For alternative flotation protocols in which (a) the cell suspension is adjusted to a density of 1.085 g/ml and the DC allowed to float up through a 1.065 g/ml density solution layered on top and (b) the cell suspension itself is adjusted to a density of 1.055 g/ml see respectively [OptiPrep™ Application Sheet C20 →→](#) and [OptiPrep™ Application Sheet C21. →→](#)

2. Preparation of cells

For the standard preparation of DCs from spleen and thymus by collagenase digestion and filtration see [OptiPrep™ Application Sheet C21. →→](#) Lung DCs are prepared by perfusion of the tissue via the pulmonary artery with phosphate-buffered saline containing heparin prior to digestion with collagenase (see ref 7. for more details).

For the preparation of DCs from bone marrow cell cultures, the cells flushed from murine femurs and tibias are cultured in the standard RPMI/10% FCS in the presence of 10 ng/ml IL-4 and granulocyte-macrophage colony-stimulating factor for 6 days prior to the harvesting of loosely adherent and non-adherent cells (see ref. 5 for more details of this cell preparation protocol).

3. Purification of dendritic cells (adapted from ref 5)

3a. Solutions required

- OptiPrep™ (shake the bottle gently before use)
- Diluent: 0.88% (w/v) NaCl, 1 mM EDTA, 0.5% bovine serum albumin (BSA), 10 mM Tricine-NaOH, pH 7.4 (see Note 1)
- Density barrier: 2.3 vol. of Solution A + 9.7 vol. of Solution B (see Note 2)
- Hank's Balanced Salt Solution (see Note 1)

Keep the following stock solutions at 4°C:
100 mM Tricine: 1.79 g per 100 ml water
100 mM EDTA (Na₂·2H₂O): 3.72g per 100 ml water

Solution B: Dissolve 0.88 g NaCl in 50 ml water; add 10 ml and 1 ml respectively of stock buffer and EDTA solutions; adjust to pH 7.4 with NaOH; dissolve 0.5 g of bovine serum albumin in the solution; check the pH and make up to 100 ml.

3b. Protocol

Carry out all operations at 4°C, making sure that all solutions and equipment are pre-cooled.

1. Harvest the cells by centrifugation at 540 g for 5 min and wash them twice in Solution D.
2. Suspend the washed cell pellet in this solution (6 ml).
3. Transfer 6 ml of Solution C to a centrifuge tube and overlay with cell suspension (see Note 3).
4. Centrifuge at 600-700 g for 5 min (see Note 5).
5. Allow the rotor to decelerate without the brake and harvest the DCs from the interface (see Figure 1 and Note 6).

4. Notes

1. The solution used to suspend the cells prior to layering on the density barrier and the solution used to dilute the OptiPrep™ may be any compatible isoosmotic medium. Hansen et al [8] used Hank's Balanced Salt Solution (HBSS) containing 5% FCS and 2 mM EDTA for both, while Cervantes-Barragan et al [6] replaced the HBSS with PBS and Luckashenak et al [7] used PBS for suspending the cells and HBSS for diluting the OptiPrep™ (both containing 0.2% BSA and 2 mM EDTA).
2. The density of this barrier is approx. 1.067 g/ml and equivalent to 11.6% (w/v) iodixanol; others have used 12% [6]. For lung DCs Hansen et al [8] employed a two-layer gradient of 4% and 16% (w/v) iodixanol.
3. The relative volumes of sample and density barrier are probably not very critical.
4. The time may be increased to 20 min. Most centrifugations are carried out at room temperature but there is no obvious reason why lower temperatures may not be used if necessary.
5. In the case of the two-layer gradient, the DCs band at the lower interface. Eaton et al [5] reported that the purity of the DCs from bone marrow cultures, as judged by flow cytometry, was approx. 85%. The harvested DC may be further purified by using the appropriate MAb-coated magnetic beads. Carrying out this initial gradient purification step allows the bead purification by negative selection to be performed more efficiently. Ref. 9 reviews DC methodology and function.

5. References

To access abstracts of refs 1-9 (file CA47) click on the double blue arrow →→

1. Barfoot, R., Denham, S., Gyure, L. A., Hall, J. G., Hobbs, Jackson, L. E. and Robertson, D. (1989) *Some properties of dendritic macrophages from peripheral lymph* Immunology, **68**, 233-239
2. McLellan, A. D., Starling, G. C. and Hart, D. N. J. (1995) *Isolation of human blood dendritic cells by discontinuous Nycodenz gradient centrifugation* J. Immunol. Meth., **184**, 81-89
3. Bonneau, M., Epardaud, M., Payot, F., Niborski, V., Thoulouze, M-I., Bernex, F., Charley, B., Riffault, S., Guilloteau, L.A. and Schwartz-Cornil, I. (2006) *Migratory monocytes and granulocytes are major lymphatic carriers of Salmonella from tissue to draining lymph node* J. Leukoc. Biol., **79**, 268-276
4. Lee, J-E., Kang, C-S., Guan, X-Y., Kim, B-T., Kim, S-H., Lee, Y-M., Moon, W-S. and Kim, D-K. (2007) *Discoidin domain receptor 2 is involved in the activation of bone marrow-derived dendritic cells caused by type I collagen* Biochem. Biophys. Res. Comm., **352**, 244-250
5. Eaton, K.A., Benson, L.H., Haeger, J. and Gray, B.M. (2006) *Role of transcription factor T-bet expression by CD4+ cells in gastritis due to Helicobacter pylori in mice* Infect. Immun., **74**, 4673-4684
6. Cervantes-Barragan, L., Züst, R., Weber, F., Spiegel, M., Lang, K.S., Akira, S., Thiel, V. and Ludewig, B. (2007) *Control of coronavirus infection through plasmacytoid dendritic cell-derived type I interferon* Blood, **109**, 1131-1137
7. Luckashenak, N.A., Ryszkiewicz, R.L., Ramsey, K.D. and Clements, J.L. (2006) *The Src homology 2 domain-containing leukocyte protein of 76-kDa adaptor links integrin ligation with p44/42 MAPK phosphorylation and podosome distribution in murine dendritic cells* J. Immunol., **177**, 5177-5185
8. Hansen, S., Lo, B., Evans, K., Neophytou, P., Holmskov, U. and Wright, J.R. (2007) *Surfactant protein D augments bacterial association but attenuates major histocompatibility complex class II presentation of bacterial antigens* Am. J. Respir. Cell. Mol. Biol., **36**, 94-102
9. Ardavin, C., Martinez del Hoyo, G., Martin, P., Anjuere, F., Arias, C. F., Marin, A. R., Ruiz, S., Parrillas, V. and Hernandez, H. (2001) *Origin and differentiation of dendritic cells* Trends Immunol., **22**, 691-700